
CAT

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Homogenising System X360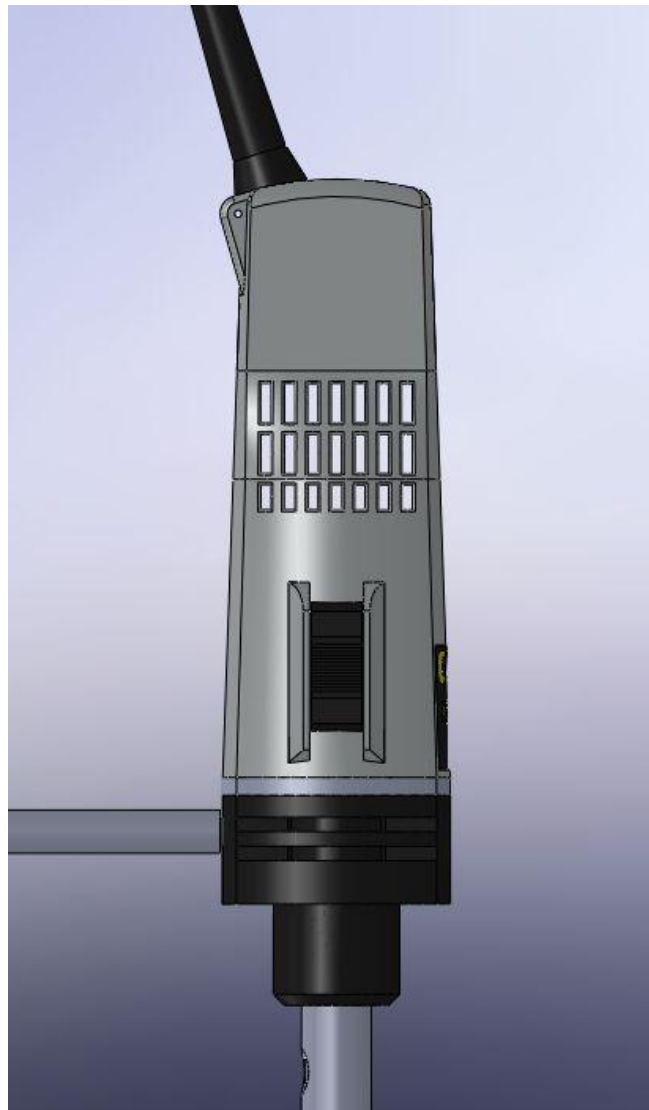

Content

1	General Information	5
2	Unpacking the Instrument	5
2.1	Scope of delivery	5
2.2	Homogenising tools and accessories (to be ordered separately):	5
3	Safety Information	7
3.1	Explanation of the safety information in the instruction manual	7
3.2	Explanation of the safety information on the device	7
3.3	General Safety Information	7
4	Set-up and Operation	8
4.1	Intended Use.....	8
4.2	Description Drive Unit X360.....	8
4.2.1	Assembly of support rod.....	9
4.2.2	Attaching the drive unit to a stand	9
4.3	Description of a homogenising tool	10
4.3.1	Inserting the homogenizing tool	10
4.4	Operation X360	11
4.4.1	Connect the unit to the mains.	11
4.5	Operation with power control unit P1500.....	12
4.5.1	Description of the power control unit P1500 for X360	12
4.5.2	Set-up of X360 with power control unit P1500	13
4.6	Exchanging the homogenizing Shafts	13
4.7	Exchanging the Generator:	13
5	Maintenance and Cleaning	14
5.1	Cleaning the drive unit X360.....	14
5.2	Cleaning the homogenising tool	14
5.3	Maintenance of drive unit X360.....	14
5.4	Maintenance homogenizing tools	14
5.5	Exchange of fuse.....	15
6	Dismantling, Transport and Storage	15
6.1	Dismantling.....	15
6.2	Transport and Storage.....	15
7	Disposal	16
8	Warranty and Liability	16
9	Technical Data.....	17

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1 General Information

The CAT homogeniser X360 is designed in accordance with Safety Class 1 and built and tested in accordance with DIN EN 61010.

According to these regulations the unit is designed to meet the requirements for safe and correct operations. To maintain the proper safety and operational functions of the instrument the user should follow the instructions and safety guidelines in this manual.

2 Unpacking the Instrument

Unpack the instrument carefully and check to see that it is not damaged. It is important that any damage incurred in transport to be recognized at the time of unpacking. Notify your carrier or forwarding agent immediately in case of such damage.

2.1 Scope of delivery

Please check that the package contains the following:

Type	Description	Part No.
X 360	Homogenizing Drive Unit, 230V	60401-00 or
X 360	Homogenizing Drive Unit, 115V	60401-01 or
	1 Support Rod	
	1 Instruction Manual	

2.2 Homogenising tools and accessories (to be ordered separately):

Tools	Description	Part No.
T 6	Shaft, 6 mm diameter, 55 mm long, complete with generator	60420*
T 6	Shaft, 6 mm diameter, 120 mm long, complete with generator	60410*
T 10	Shaft, 10 mm diameter complete with generator	60421*
T 17	Shaft, 10 mm diameter, complete with 17 mm generator	60426*
T 20	Shaft, 20 mm diameter complete with generator	60422*
G 20	Shaft, 20 mm diameter complete with generator	60423*
Pedestal stand incl. rod	platform overall dimensions: 330x200 mm, rod diameter: 16mm, length:690mm	60491-00
Cross over clamp	Casted alloy clamp to attach the unit to a stand	60492-00
Cross over clamp	Steel clamp to attach the unit to a stand	60492-50
Universal key	Universal rotor wrench for Shafts 6 - 30 mm	60470-00
Socket wrench	for all shafts	60471-00
Rotor wrench	for shaft T 10	60472-00

**Important Note:**

Read this instruction manual carefully before operating the instrument. Should there be any additional questions, after reading these instructions, concerning the set-up, operation or warranty, please contact either your distributor, or the manufacturer at the following address

CAT

Ingenieurbüro CAT
M. Zipperer GmbH
Etzenbach 16
D-79219 Staufen
+49-7636-7803-0

**Important Note:**

Please store the instruction manual in a place easily accessible to every user.

**Important Note:**

Make sure that the unit is to be operated under the conditions corresponding to that as described under the section titled "Technical Data"

A EURO AC plug (DIN 49441 CEE 7/VII10/ 16 A 250 V) is standard on all instruments. For North America instruments feature standard US plugs (NEMA Pub.No.WDI1961 ASA C 73.1. 1961 page 8 15A 125V).

When operating the instruments in countries with different AC plug systems use an approved adapter or have a qualified electrician replace the AC plug with an approved model suitable for the country of operation.

The instrument is earthed as supplied. When replacing the original AC plug, ensure that the earth conductor is connected to the new plug!

3 Safety Information

3.1 Explanation of the safety information in the instruction manual

This symbol emphasizes the following notices, which have to be strictly observed by the user! Disregarding of warnings may result in impairment of serviceability as well as impairment of the user.

This symbol emphasizes prohibitions, which have to be strictly observed. Any disregard of the prohibitions may result in severe impairment of serviceability as well as severe impairment of the user.

This symbol emphasizes notices, which have to be observed by the user to assure a safe operation of the unit.

CAUTION:
Hazard of fire or explosion!

3.2 Explanation of the safety information on the device

CAUTION:
This symbol indicates to read the instruction manual carefully prior to operation of the instrument. Please mark points which require special attention in your field of application so they are not overlooked. Disregarding of warnings may result in impairment of serviceability as well as impairment of the user.

3.3 General Safety Information

Attention:
Please comply with all safety and accident-prevention regulations applicable to laboratory work.

Attention:
Do not use the unit with flammable or explosive substances. The instrument is not explosion-proof.

Attention:
Only instructed users may operate the instrument.

Important Note:
Make sure that all safety information of the instrument respectively on the device itself are clearly visible during operation

Attention:
If noise level exceeds 85 dB (A) at the work station operator must use ear plug and/or ear mufflers.

Attention:
Do not open the instrument. Repairs are only to be carried out by trained service technicians.



Ensure that the unit is firmly attached to a stand.

4 Set-up and Operation


4.1 Intended Use

The X360 is a homogenizing system which in connection with a homogenizing tool produces emulsions and dispersions. For safe operation the unit has to be fixed to a stand.

4.2 Description Drive Unit X360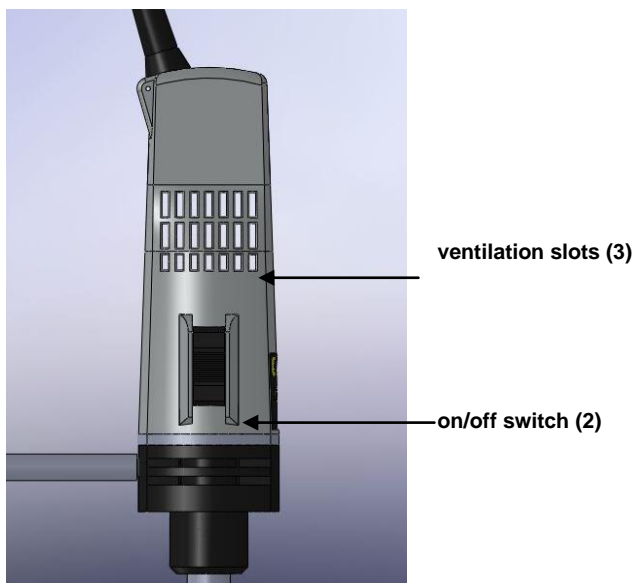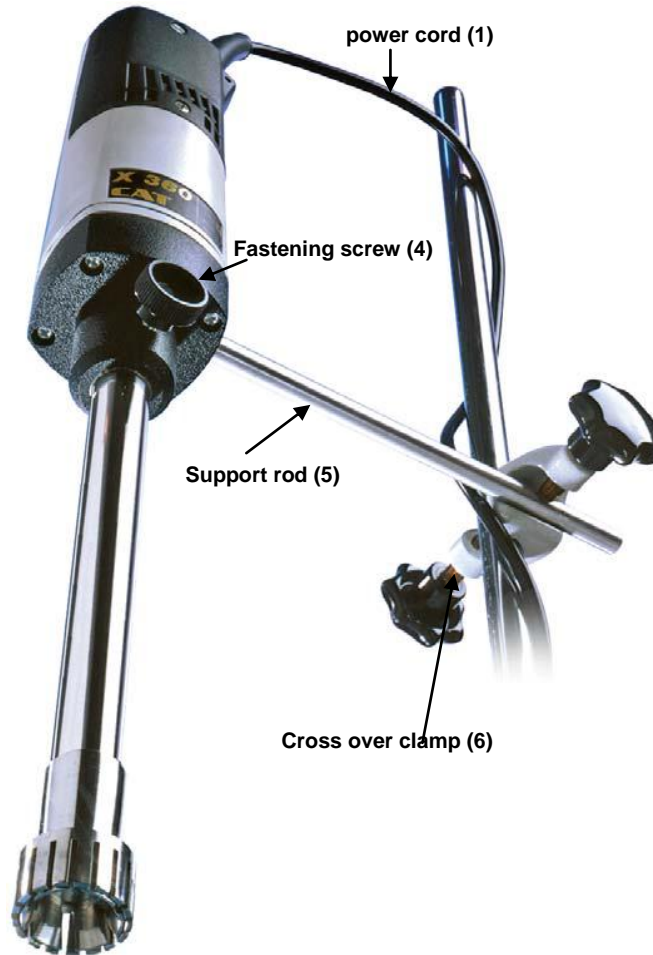

4.2.1 Assembly of support rod

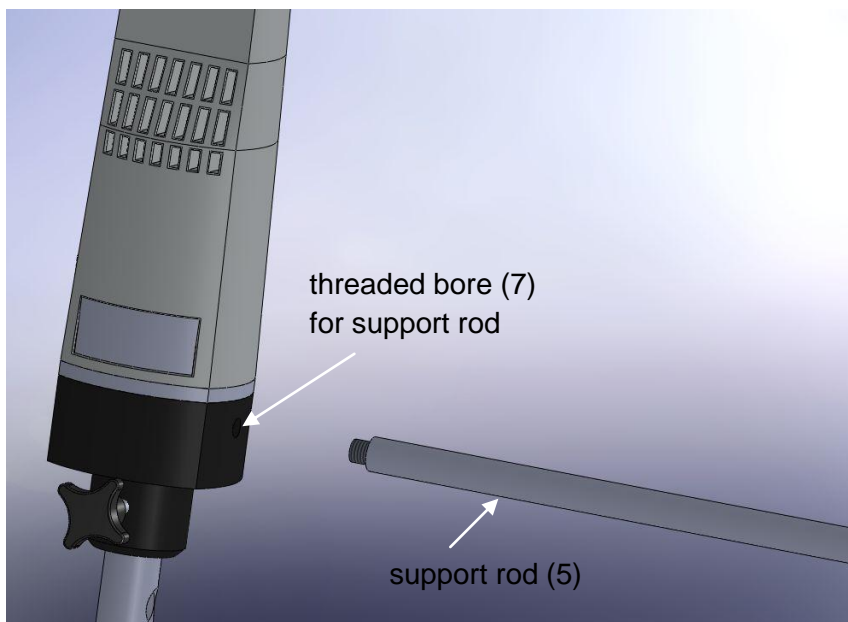
- Screw the support rod (5) into the threaded bore (7) at the back of the unit.
- Insert a screw driver into the bore at the end of the support rod (5) and firmly tighten the support rod by turning the screw driver to the right.

4.2.2 Attaching the drive unit to a stand


To ensure safe operation mount the drive unit to a stand using a cross over clamp (6). Vibrations may loosen the clamp screws (8). Please check periodically whether all fastening screws are securely tight.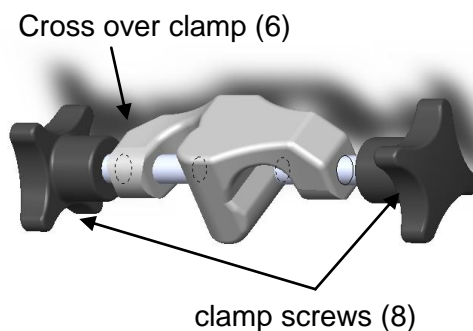

- Attach cross over clamp (6) to a stand rod.
- Now attach the drive unit to the stand by inserting the support rod (5) into the free opening of the cross over clamp (6).
- Tighten screws (8) of the cross over clamp (6).

4.3 Description of a homogenising tool

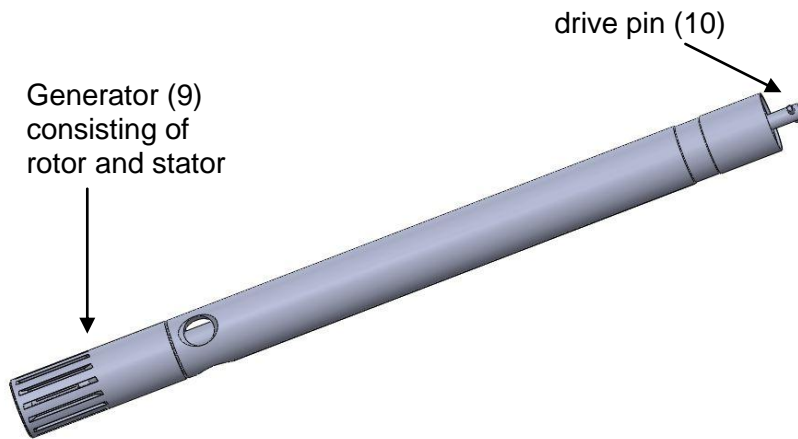
4.3.1 Inserting the homogenizing tool

The Generators (9) have sharp edges. Handle with care. Risk of injuries!

The homogenizing tool comes always assembled. The connecting adapter (drive pin (10)) is located on the top of the Shaft and is used to attach the Shaft to the drive motor. When inserting the homogenising tool, make sure that the coupling fits well. This can be easily detected by the "click" that is heard when the coupling Shaft has reached the end of the coupling slot of the motor. Securely fasten the fastening screw (4).

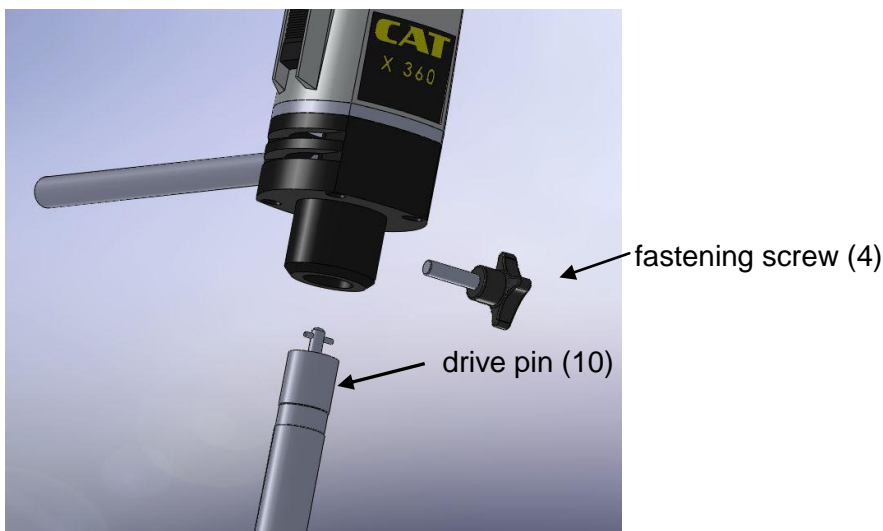
Never run homogenising tools dry as the bearings and gaskets will be damaged if the Generator is not cooled by the medium. The distance between the homogenising tool and the vessel bottom should not be less than 10mm. The filling of the medium may not be less than about 55 mm. To avoid a vortex it is necessary to insert the dispersing Shaft out of the center of the vessel. If these conditions are met the unit is ready for operation.

4.4 Operation X360

Attention:

When connecting the instrument to an AC power outlet, ensure that your local supply voltage matches that indicated on the instrument's rating plate.

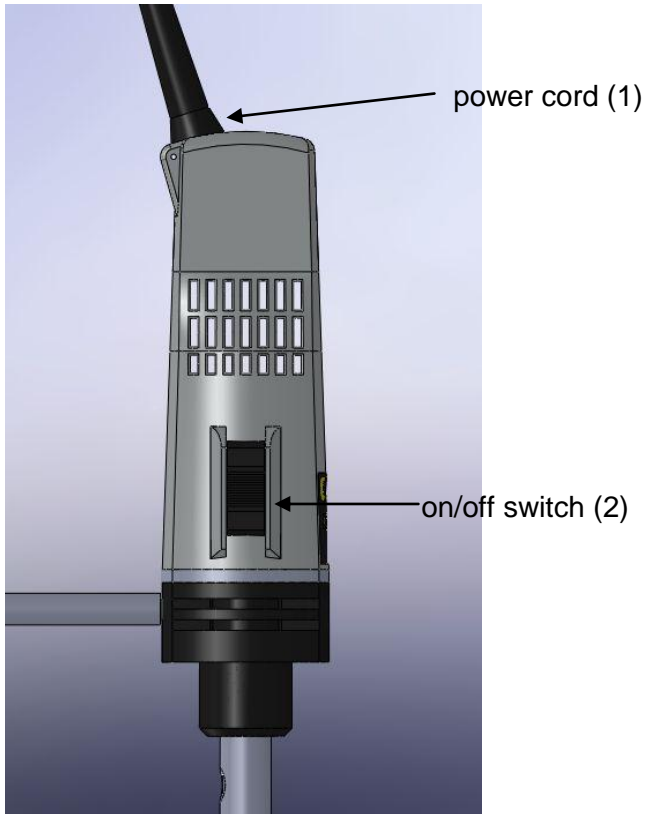4.4.1 Connect the unit to the mains.

The drive unit is turned on and off by means of the on/off switch (2) at the front of the motor housing.

position I: motor runs

position 0: motor off

Caution:

The on/off switch (2) is only single-pole and is thus not suitable for safely isolating the unit from the mains supply!

Before starting the unit make sure that

- the Generator (9) of the homogenising tool is inserted in the fluid. To avoid a vortex it is necessary to insert the dispersing Shaft out of the centre of the vessel.
- the vessel is protected against rotating, shaking or moving
- make sure that the air ventilation slots (3) are always open, as this is necessary for cooling purposes

4.5 Operation with power control unit P1500

4.5.1 Description of the power control unit P1500 for X360

The power control unit is used for continuous speed control of the X360.

 The below stated speed rates are only approximate values.

Approximate Speed Rates

Pos.	1	2	3	4	5	6	7	8	9	10
rpm	10800	13500	18000	21500	25800	30500	32800	36500	40000	42500

4.5.2 Set-up of X360 with power control unit P1500

Plug the power cable of X360 (81) into the socket (13) of the power control unit P1500

Now plug in the power cable (12) of the P1500 into a power source.

Switch on the X360

Start the unit by switching the potentiometer (14) to position "1"

The power control unit P1500 is now operating

With the potentiometer (14) the voltage can be controlled.

Always start with the lowest speed (Position "1") and then increase until the desired speed setting has been reached.

4.6 Exchanging the homogenizing Shafts

- Prior to exchange the homogenizing shaft disconnect the unit from the mains supply by unplugging the unit.
- Loosen the fastening screw (4) and firmly pull out the Shaft by hand. It will disconnect easily.
- When inserting another Shaft, make sure that the coupling fits well. This can be easily detected by the "click" that is heard when the coupling Shaft has reached the end of the coupling slot of the motor.
- Securely fasten the fastening screw (4).

4.7 Exchanging the Generator:

The Generators have sharp edges. Handle with care. Risk of injuries!

To exchange the Generator (rotor and stator), always take the Shaft out of the motor drive unit as described under 4.6.

For undoing the rotor, use the two special keys (to be ordered separately). The socket wrench is used to counter-hold the Shaft, and the rotor wrench to undo the rotor.

To exchange the stator, remember that the fastening thread is **counter-clockwise**.

You will find further detailed descriptions in the instruction manual which comes with each homogenising tool.

Important Note

Always unscrew the rotor first before unscrewing the stator.

5 Maintenance and Cleaning

5.1 Cleaning the drive unit X360

To clean the drive unit use only water with a detergent that contains tensides or use isopropyl alcohol for stubborn soiling.

5.2 Cleaning the homogenising tool

The Generators (9) have sharp edges. Handle with care. Risk of injuries!

To avoid clogging of the shaft, clean shaft, generator and in case of G-shaft the sealing parts after each use. This is done by operating it in a solvent which dissolves substance residues and is not harmful to the gasket. This is usually sufficient to clean the Generator.

Chemical sterilization may be also a method. General-purpose disinfectants such as formalin, alcohol, etc. may be used. It is important to remove disinfectant residues with sterilized water.

Important Note:

Make sure that the bearings, O-rings and gaskets are resistant to solvents.

Sterilization with moist heat:

This denotes the use of a steam jet pressurized to 2 bars at 120°C.

5.3 Maintenance of drive unit X360

Attention:

Do not open the instrument. Repairs are only to be carried out by trained service technicians.

5.4 Maintenance homogenizing tools

The Generators have sharp edges. Handle with care. Risk of injuries!

The gaskets in the homogenizing tools must be constantly monitored. In the event of leakage the suction effect of the rotating Shaft can cause the medium to penetrate as far as the drive unit. If liquid emerges from the side hole at the top of the Shaft tube stop work immediately and check the gaskets. The function of the Generators depends on the condition of the sharp edges on the rotor and stator. These edges may be blunted very quickly in abrasive media, reducing the effectiveness of homogenizing.

5.5 Exchange of fuse

- Unplug power cord (11)
- Unplug power cord (2) of X360 from socket (13) of P1500
- Open fuse (12) holder with a coin or screwdriver
- Replace fuse

Only replace fuses which are equivalent to the original fuse.

6 Dismantling, Transport and Storage

6.1 Dismantling

1. Switch the unit off.
2. Disconnect it from the mains supply.
3. Remove any glass beakers and any other equipment around the drive unit.
4. Remove the homogenizing tool by loosening the fastening screw (4).
5. Loosen the special clamp and remove the clamp from the support rod.
6. Unscrew the support rod.
7. Now you may remove the instrument from the working area.

6.2 Transport and Storage

Prior to transport:

Switch the instrument off and proceed with dismantling as described under “Dismantling”.

Place the instrument and its parts in its original packaging or another suitable container to protect it during transport. Close the packaging with adhesive tape.

Store the instrument in a dry environment. Please observe the specified conditions of the ambient (temperature and humidity).

Do not subject the instrument to mechanical shocks or vibration during transporting it.

In case you do not use the original packaging please mark the box with the following notes:

- Glass symbol (handle with care, fragile)
- Umbrella (keep dry)
- Content (list of content)
- Storage ambient:
 - Max. ambient temperature : RT to +40°C
 - Max. humidity: 80%

7 Disposal

Please dispose of used instruments and defective components at your local recycling collection point. Prior to disposal, sort according to materials: Metal, glass, plastic, etc. Also be sure to dispose of the packing material in an environmental-friendly manner.

8 Warranty and Liability

The manufacturer agrees to either repair, or replace, at the manufacturer's discretion, any defects in materials or workmanship which develop within 24 months of the delivery of this product to the original user. In the event of replacement, the replacement unit will be guaranteed for the remainder of the original twenty-four (24) months period or ninety (90) days, whichever is longer.

If this product should require service, contact your local distributor or manufacturer for necessary instructions.

This guarantee will not apply if the defect or malfunction was caused by accident, neglect, unreasonable use or fitness for a particular purpose, which extend beyond the description and period set forth herein.

The manufacturer's sole obligation under this guarantee is limited to the repair or replacement of a defective product and the manufacturer shall not, in any event, be liable for any incidental or consequential damages of any kind, resulting from use or possession of the product.

Attention:

The user has to determine, if the instrument is suitable for his specific application. If there are any further queries, contact your local dealer or the manufacturer direct.

9 Technical Data

9.1 Technical Data X360

Type: X360	Specifications
Motor Power intake	350 Watt
Idle speed	Fixed 45000 rpm
With P1500	10.800 – 42500 rpm
Voltage	230V/50 Hz or 110V/60 Hz
Case dimensions	290x63x62 mm, LxWxD
Weight	1.3 kg

9.2 Technical Data P1500

Type: P1500	Specifications
Power Supply	230 V AC 50/60 Hz, 1500 W
Output	25 – 228 V AC stepless and loss-free
Electronics	Phase-angle control
Connection	Cable, 120 cm long with grounded plug
Load	Earthed socket (Schuko)
Case dimensions	150x80x55 mm, LxWxH
Weight	0.5 kg
Fuse	6.3A, FF 250V, 5x20mm
X360 and P1500	
Permissible ambient temperature	5 - 40 °C
Permissible humidity	80 % RH

10 Repairs

When returning instruments for repair that have come into contact with hazardous substances, please:

Fill in attached "Repair Return Form"

Provide precise information on the relevant medium

Take protective measures to ensure the safety of our receiving and maintenance personnel

Mark the package as appropriate for hazardous materials.

10.1 Repair Return Form

CONTACT/USER INFORMATION

Contact:	Phone No.
Fax No.	Email:
Billing:	Shipping:
Company	Company
Address	Address

INSTRUMENT INFORMATION

Model	Serial No.	
Please describe all problems/malfunctions		
Operating Conditions (please fill in if applicable)		
Ambient Temp.	Humidity	Speed
Load	Volume	Viscosity
Temperature in °C	Sample Temperature	Operating Time
Sample Description*		

*NOTE: If the instrument was exposed to hazardous material, it must be decontaminated BEFORE returning it to CAT and an MSDS for hazardous material must be included with the instrument.

RETURN SHIPPING

UPS <input type="checkbox"/>	Air Parcel Post <input type="checkbox"/>	Collect** <input type="checkbox"/>	Other** <input type="checkbox"/>
------------------------------	--	------------------------------------	----------------------------------

**Your account number is required for UPS collect respectively the address and contact of your preferred forwarder if you choose any other transport means.

PACKAGING INSTRUCTIONS TO RETURN A CAT INSTRUMENT FOR REPAIR

- ✓ Remove all accessories (e.g. homogniser tools, stirring paddles) from the instrument
- ✓ Clean excess testing material off the instrument/accessory
- ✓ Include MSDS sheets for all hazardous materials used with this instrument
- ✓ Pack the instrument in its original box. If the box is not available, take care to wrap the instrument and accessories with enough material to support them.
- ✓ DO NOT send pedestal stand unless there is a problem with the upright rod, clamp or base. If there is a problem with the stand remove the upright rod from the base and individually wrap each item to avoid contact with the instrument. (Applicable for overhead stirrers and homogenisers.)
- ✓ Pack the instrument and related items in a strong box for shipping. Mark the outside of the box with handling instructions.

Example: "Handle with care" or "Fragile- Delicate Instrument"

Ingenieurbüro
M. Zipperer GmbH

Etzenbach 16,D-79219 Staufen
Tel.: +49 (0) 7636 78030
Fax: +49 (0) 7636 780345
E-Mail: info@cat-ing.de
Internet: <http://www.cat-ing.de>

CE Declaration of conformity

We declare under our sole responsibility that this product (see label) corresponds to the regulations and conforms to the standards or standardized documents

2006/95/EC

DIN EN 61010-1; VDE 0411-1:2002-08;Amendment 1; VDE 0411-1 Amendment 1:2002-11; Amendment 2; VDE 0411-1 Amendment 2:2004-01
DIN EN 61010-2-051; VDE 0411-2-051:2004-07

2004/108/EC

DIN EN 61326-1; VDE 0843-20-1:2006-10
DIN EN 61000-6-2; VDE 0839-6-2:2006-03
DIN EN 61000-6-4; VDE 0839-6-4:2007-09
DIN EN 61000-4-2; VDE 0847-4-2:2009-12
DIN EN 61000-4-3; VDE 0847-4-3:2008-06
DIN EN 61000-4-4; VDE 0847-4-4:2005-07
DIN EN 61000-4-5; VDE 0847-4-5:2007-06
DIN EN 61000-4-6; VDE 0847-4-6:2009-12
DIN EN 55011; VDE 0875-11:2010-05
DIN EN 55014-1; VDE 0875-14-1:2010-02
DIN EN 55014-2; VDE 0875-14-2:2009-06

2006/42/EC, 93/68/EEC

DIN EN ISO 12100-1:2004-04
DIN EN ISO 12100-2:2004-04

In the case of a modification of the unit which has not been agreed on with us, this declaration becomes null and void and the warranty expires.

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