**CAT**

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## *Instruction Manual*

---

### *Homogenising System UNIDRIVE X1000 and X1000D*

## Content

1	General Information .....	5
2	Unpacking the Instrument .....	5
2.1	Scope of Delivery.....	5
2.2	Homogenising Tools and Accessories (to be ordered separately): .....	5
3	Safety Information .....	7
3.1	Explanation of the Safety Information in the Instruction Manual .....	7
3.2	Explanation of the Safety Information on the Device .....	8
3.3	General Safety Information .....	8
4	Set-up and Operation .....	8
4.1	Intended Use.....	8
4.2	Description Drive Unit UNIDRIVE X1000/X1000D.....	9
4.3	Set-up .....	10
4.3.1	Assembly of Support Rod.....	10
4.3.2	Attaching the Drive Unit to a Stand .....	11
4.4	Description of a Homogenising Tool .....	11
4.4.1	Inserting the Homogenizing Tool .....	12
4.5	Operation UNIDRIVE X1000/X1000D .....	13
4.5.1	Connect the Unit to the Mains. ....	13
4.5.2	Description of the Speed Scale X1000 and X1000D: .....	14
4.6	Exchanging the Homogenizing Shafts .....	14
4.6.1	Exchanging the Generator:.....	14
5	Assembling the DK 30 Flow-Through-Chamber.....	15
5.1	Description of DK 30 Flow-Through-Chamber .....	15
5.2	Mounting the Flow-Through-Chamber to the Drive Unit UNIDRIVE X1000/X1000D ...	16
6	Operation of Flow Through-Chamber.....	17
6.1	Flow-Through Mode .....	17
6.2	Circulation Mode .....	17
7	Analytical Mill Attachment AX 60.....	17
7.1	Description of the Analytical Mill Attachment AX 60 .....	17
8	Mounting of the Analytical Mill AX 60 .....	18
8.1	Exchange of the cutting knife .....	19
9	Maintenance and Cleaning .....	19
9.1	Cleaning the Drive Unit UNIDRIVE X1000/X1000D .....	19
9.2	Cleaning the Homogenising Tool.....	19

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## 1 General Information

The CAT homogeniser Unidrive X1000/Unidrive X1000D is designed in accordance with Safety Class 1 and built and tested in accordance with DIN EN 61010.

According to these regulations the unit is designed to meet the requirements for safe and correct operations. To maintain the proper safety and operational functions of the instrument the user should follow the instructions and safety guidelines in this manual.

## 2 Unpacking the Instrument

Unpack the instrument carefully and check to see that it is not damaged. It is important that any damage incurred in transport to be recognized at the time of unpacking. Notify your carrier or forwarding agent immediately in case of such damage.

### 2.1 Scope of Delivery

Please check that the package contains the following:

Type	Description	Part No.
X 1000	Homogenizing Drive Unit, 230V	60402-50 or
X 1000	Homogenizing Drive Unit, 115V	60402-51 or
X 1000D	Homogenizing Drive Unit, 230V	60406-50 or
X 1000D	Homogenizing Drive Unit, 115V	60406-51 or
	1 Support Rod	
	1 Instruction Manual	

### 2.2 Homogenising Tools and Accessories (to be ordered separately):

Tools	Description	Part No.
T 6	Shaft, 6 mm diameter, 55 mm long, complete with generator	60420*
T 6	Shaft, 6 mm diameter, 120 mm long, complete with generator	60410*
T 10	Shaft, 10 mm diameter complete with generator	60421*
T 17	Shaft, 10 mm diameter, complete with 17 mm generator	60426*
T 20	Shaft, 20 mm diameter complete with generator	60422*
G 20	Shaft, 20 mm diameter complete with generator	60423*
T 30	Shaft, 20 mm diameter complete with 30 mm generator	60424*
G 30	Shaft, 20 mm diameter complete with 30 mm generator	60425*
DK 30	Flow-through chamber for shaft G20 or G30	60428-00
AX60	Analytical mill	60409-00

<b>Tools</b>	<b>Description</b>	<b>Part No.</b>
<b>Spare knife</b>	Knife for AX60 made of stainless steel	60478-00
<b>Spare knife</b>	Knife for AX60 made of stainless steel with carborundum insert	60479-00
<b>Pedestal stand incl. rod</b>	platform overall dimensions: 330x200 mm, rod diameter: 16mm, length:690mm	60491-00
<b>Cross over clamp</b>	Casted alloy clamp to attach the unit to a stand	60492-00
<b>Cross over clamp</b>	Steel clamp to attach the unit to a stand	60492-50
<b>Universal key</b>	Universal rotor wrench for Shafts 6 - 30 mm	60470-00
<b>Socket wrench</b>	for all shafts	60471-00
<b>Rotor wrench</b>	for shaft T 10	60472-00

**Important Note:**

Read this instruction manual carefully before operating the instrument. Should there be any additional questions, after reading these instructions, concerning the set-up, operation or warranty, please contact either your distributor, or the manufacturer at the following address

# CAT

Ingenieurbüro CAT  
M. Zipperer GmbH  
Etzenbach 16  
D-79219 Staufen  
+49-7636-7803-0

**Important Note:**

Please store the instruction manual in a place easily accessible to every user.

**Important Note:**

Make sure that the unit is to be operated under the conditions corresponding to that as described under the section titled "Technical Data"

A EURO AC plug (DIN 49441 CEE 7/VII10/ 16 A 250 V) is standard on all instruments. For North America instruments feature standard US plugs (NEMA Pub.No. WDI1961 ASA C 73.1. 1961 page 8 15A 125V).

When operating the instruments in countries with different AC plug systems use an approved adapter or have a qualified electrician replace the AC plug with an approved model suitable for the country of operation.

The instrument is earthed as supplied. When replacing the original AC plug, ensure that the earth conductor is connected to the new plug!

## 3 Safety Information

### 3.1 Explanation of the Safety Information in the Instruction Manual

This symbol emphasizes the following notices, which have to be strictly observed by the user! Disregarding of warnings may result in impairment of serviceability as well as impairment of the user.

This symbol emphasizes prohibitions, which have to be strictly observed. Any disregard of the prohibitions may result in severe impairment of serviceability as well as severe impairment of the user.

This symbol emphasizes notices, which have to be observed by the user to assure a safe operation of the unit.

**CAUTION:**

Hazard of fire or explosion!

### 3.2 Explanation of the Safety Information on the Device

**CAUTION:**

*This symbol indicates to read the instruction manual carefully prior to operation of the instrument. Please mark points which require special attention in your field of application so they are not overlooked. Disregarding of warnings may result in impairment of serviceability as well as impairment of the user.*

### 3.3 General Safety Information

**Attention:**

*Please comply with all safety and accident-prevention regulations applicable to laboratory work.*

**Attention:**

*Do not use the unit with flammable or explosive substances. The instrument is not explosion-proof.*

**Attention:**

*Only instructed users may operate the instrument.*

**Important Note:**

*Make sure that all safety information of the instrument respectively on the device itself are clearly visible during operation*

**Attention:**

*If noise level exceeds 85 dB (A) at the work station operator must use ear plug and/or ear muffers.*

**Attention:**

*Do not open the instrument. Repairs are only to be carried out by trained service technicians.*

*Ensure that the unit is firmly attached to a stand.*

*Wear ear protection should the noise level exceed 85 dB(A).*

*Wear eye protection.*

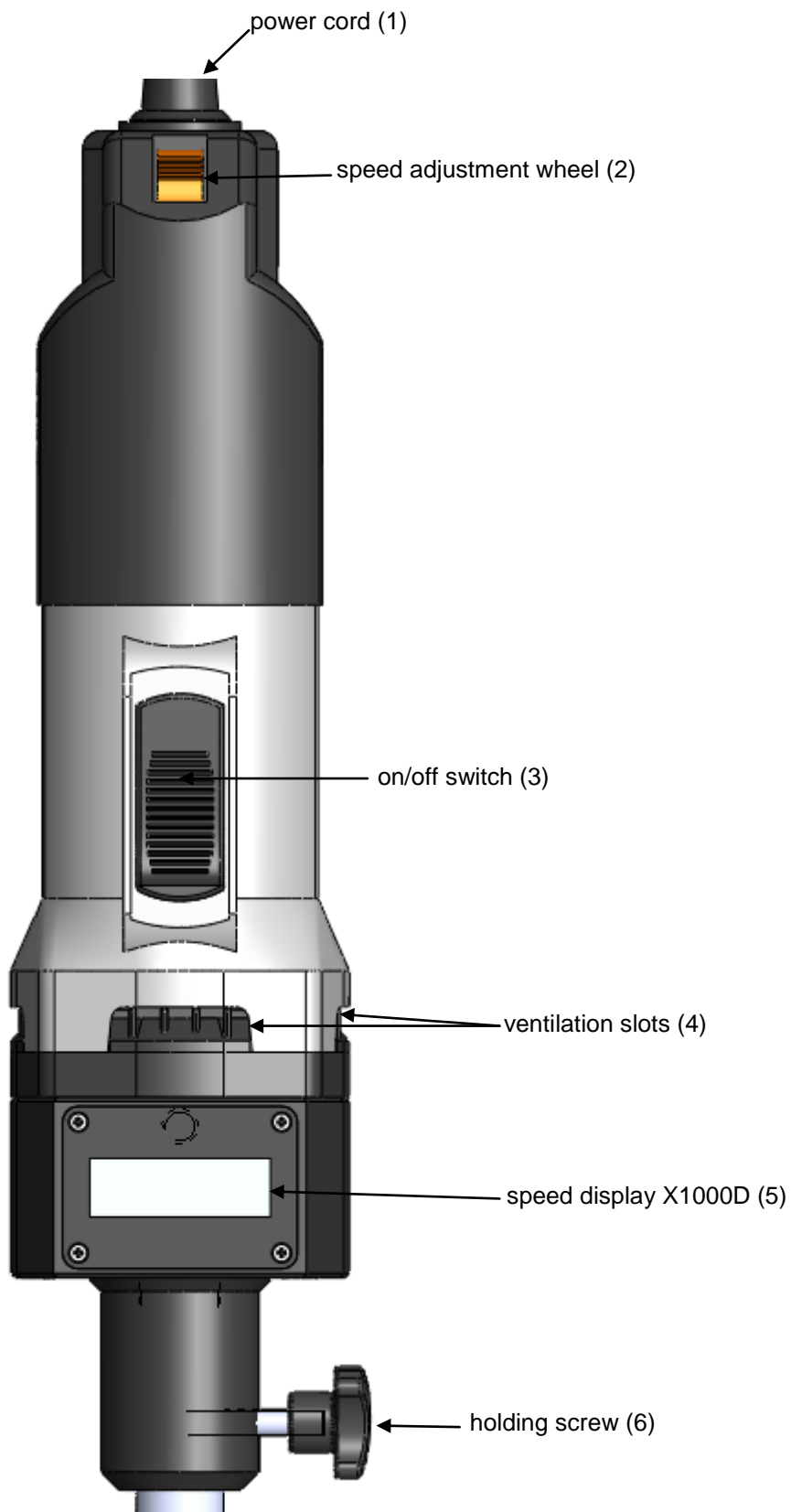
## 4 Set-up and Operation

### 4.1 Intended Use

The UNIDRIVE X1000/X1000D is a homogenizing system which in connection with a homogenizing tool produces emulsions and dispersions. For safe operation the unit has to be fixed to a stand.


## 4.2 Description Drive Unit UNIDRIVE X1000/X1000D

## 4.3 Set-up
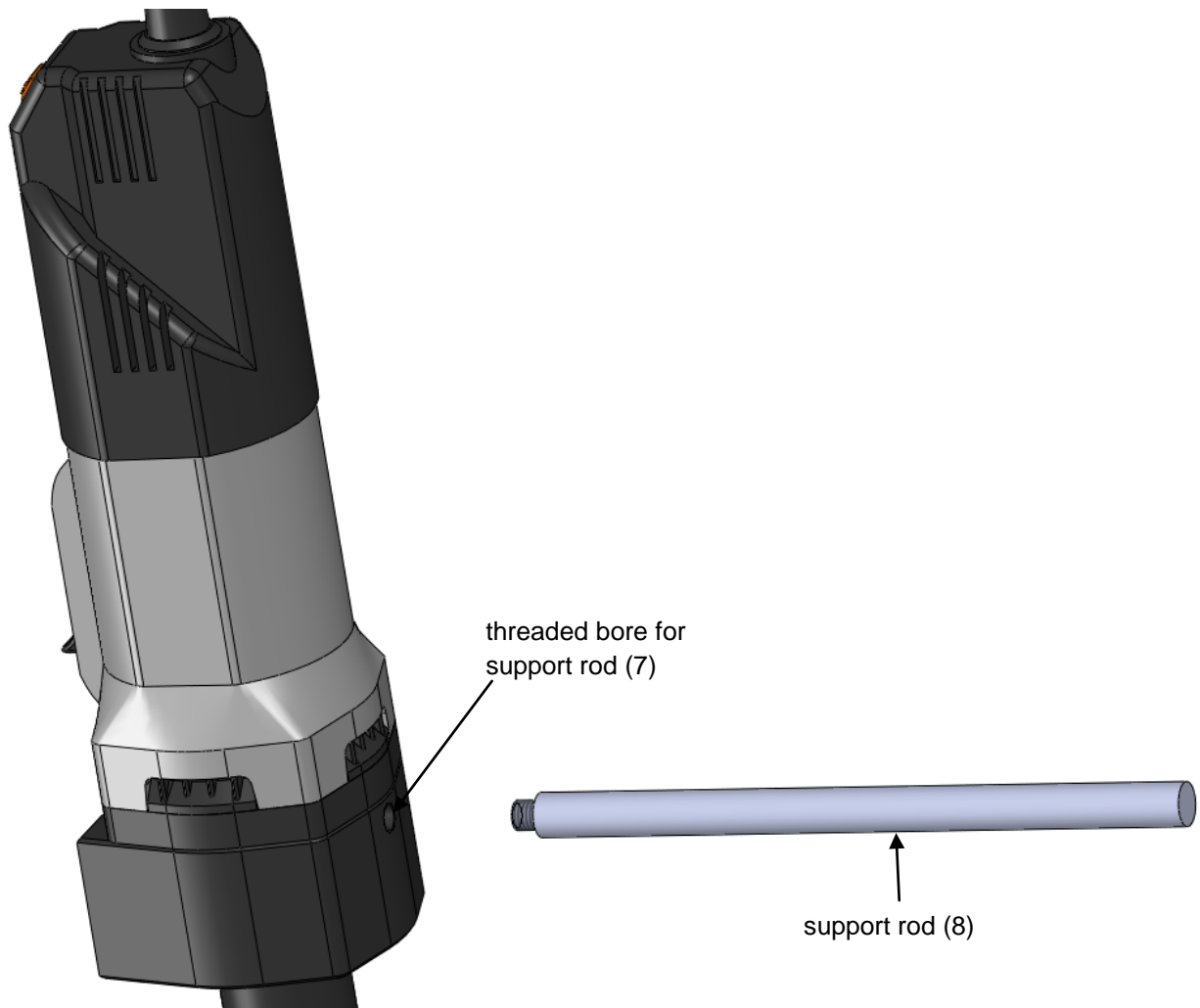
### 4.3.1 Assembly of Support Rod


- Screw the support rod (8) into the threaded bore (7) at the back of the unit.
- Insert a screw driver into the bore at the end of the support rod (8) and firmly tighten the support rod by turning the screw driver to the right.

### 4.3.2 Attaching the Drive Unit to a Stand

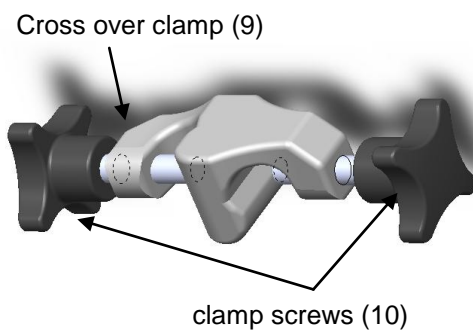
To ensure safe operation mount the drive unit to a stand using a cross over clamp (9). Vibrations may loosen the clamp screws (10). Please check periodically whether all holding screws are securely tight.

- Attach cross over clamp (9) to a stand rod.
- Now attach the drive unit to the stand by inserting the support rod (8) into the free opening of the cross over clamp (9).
- Tighten screws (10) of the cross over clamp (9).

### 4.4 Description of a Homogenising Tool

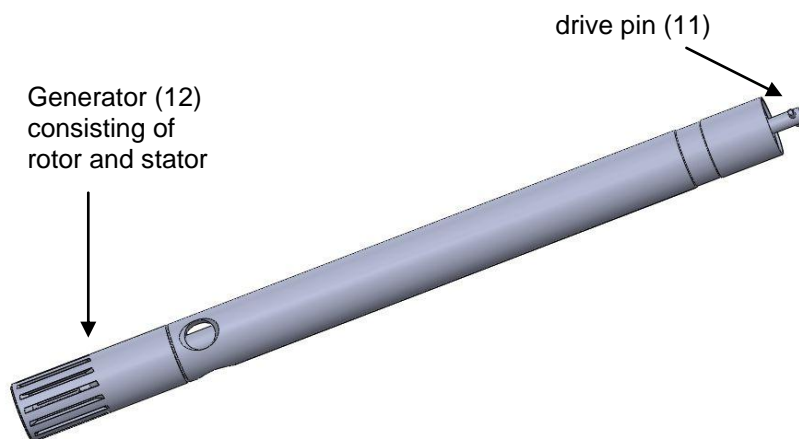
#### 4.4.1 Inserting the Homogenizing Tool

*The Generators have sharp edges. Handle with care. Risk of injuries!*

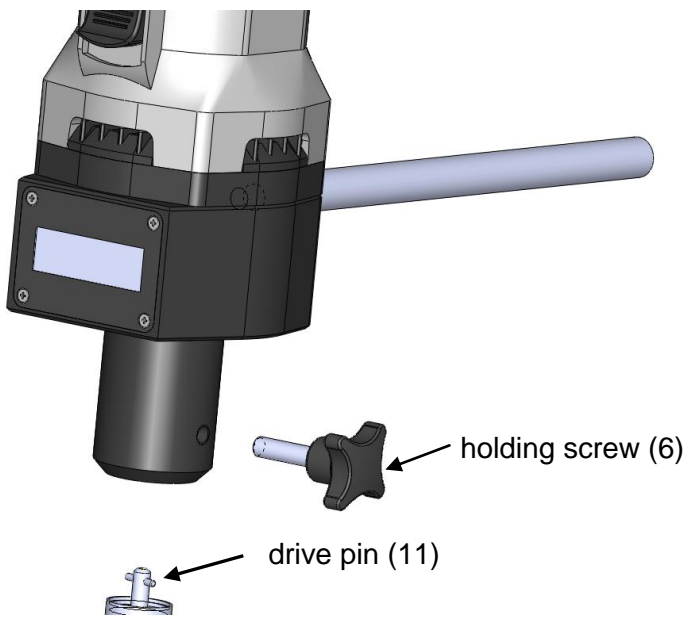The homogenizing tool comes always assembled. The connecting adapter (drive pin (11)) is located on the top of the Shaft and is used to attach the Shaft to the drive motor. When inserting the homogenising tool, make sure that the coupling fits well. This can be easily detected by the "click" that is heard when the coupling Shaft has reached the end of the coupling slot of the motor. Securely fasten the holding screw (6).


***Never run homogenising tools dry as the bearings and gaskets will be damaged if the Generator is not cooled by the medium. The distance between the homogenising tool and the vessel bottom should not be less than 10mm. The filling of the medium may not be less than about 55 mm. To avoid a vortex it is necessary to insert the dispersing Shaft out of the centre of the vessel. If these conditions are met the unit is ready for operation.***

## 4.5 Operation UNIDRIVE X1000/X1000D

### **Attention:**

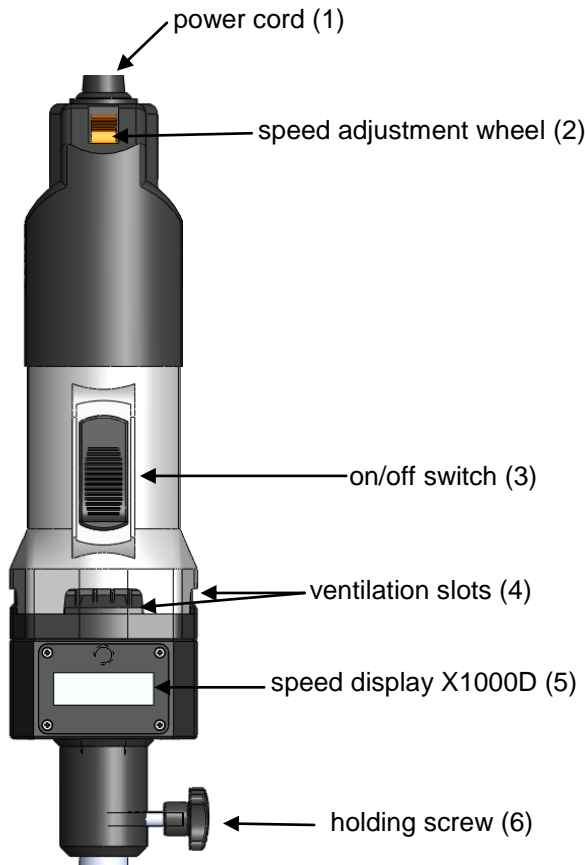*When connecting the instrument to an AC power outlet, ensure that your local supply voltage matches that indicated on the instrument's rating plate.*

### 4.5.1 Connect the Unit to the Mains.

The drive unit is turned on and off by means of the on/off switch (3) at the front of the motor housing.

position I: motor runs  
position 0: motor off

Set the speed with the speed adjustment wheel (2).

### **Caution:**

*The on/off switch (3) is only single-pole and is thus not suitable for safely isolating the unit from the mains supply!*

Before starting the unit make sure that

- the speed is set to the lowest speed
- the Generator (12) of the homogenising tool is inserted in the fluid. To avoid a vortex it is necessary to insert the dispersing Shaft out of the centre of the vessel.
- the vessel is protected against rotating, shaking or moving
- make sure that the air ventilation slots (4) are always open, as this is necessary for cooling purposes

Start the unit always at the lowest speed and then gradually increase the speed to the desired level.

#### 4.5.2 Description of the Speed Scale X1000 and X1000D:

1	=	4000 rpm
2	=	8500 rpm
3	=	15000 rpm
4	=	20000 rpm
5	=	25000 rpm
6	=	33000 rpm

The set speed of UNIDRIVE X1000D is shown the speed display (5).

#### 4.6 Exchanging the Homogenizing Shafts

- Prior to exchange the homogenizing shaft disconnect the unit from the mains supply by unplugging the unit.
- Loosen the holding screw (6) and firmly pull out the Shaft by hand. It will disconnect easily.
- When inserting another Shaft, make sure that the coupling fits well. This can be easily detected by the "click" that is heard when the coupling Shaft has reached the end of the coupling slot of the motor.
- Securely fasten the holding screw (6).

##### 4.6.1 Exchanging the Generator:

*The Generators have sharp edges. Handle with care. Risk of injuries!*

To exchange the Generator (rotor and stator), always take the Shaft out of the motor drive unit as described under 4.6.

For undoing the rotor, use the two special keys (to be ordered separately). The socket wrench is used to counter-hold the Shaft, and the rotor wrench to undo the rotor.

To exchange the stator, remember that the fastening thread is **counter-clockwise**.

You will find further detailed descriptions in the instruction manual which comes with each homogenising tool.

#### **Important Note**

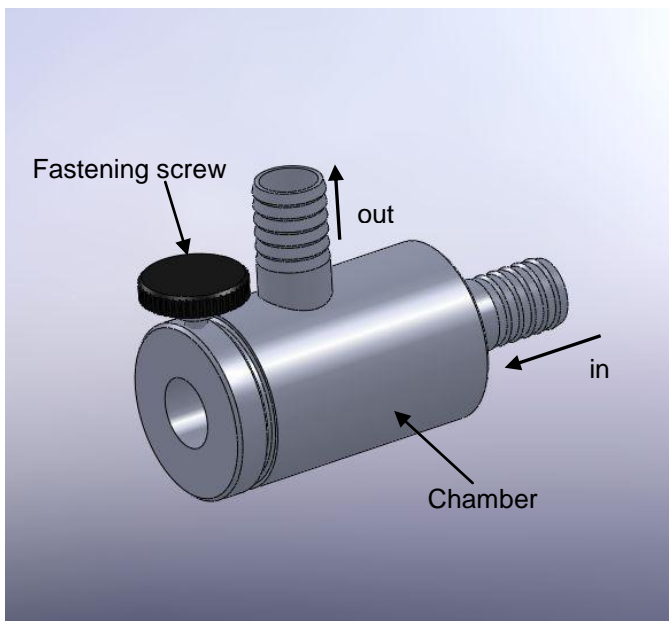
*Always unscrew the rotor first before unscrewing the stator.*

## 5 Assembling the DK 30 Flow-Through-Chamber


### 5.1 Description of DK 30 Flow-Through-Chamber

The Flow-Through-Chamber DK 30 plus the G20 or G30 dispersing Shaft makes it possible to disperse on through-flow basis.

The results are the same as those attained with the G20 or G30 on a batch basis.

## 5.2 Mounting the Flow-Through-Chamber to the Drive Unit UNIDRIVE X1000/X1000D

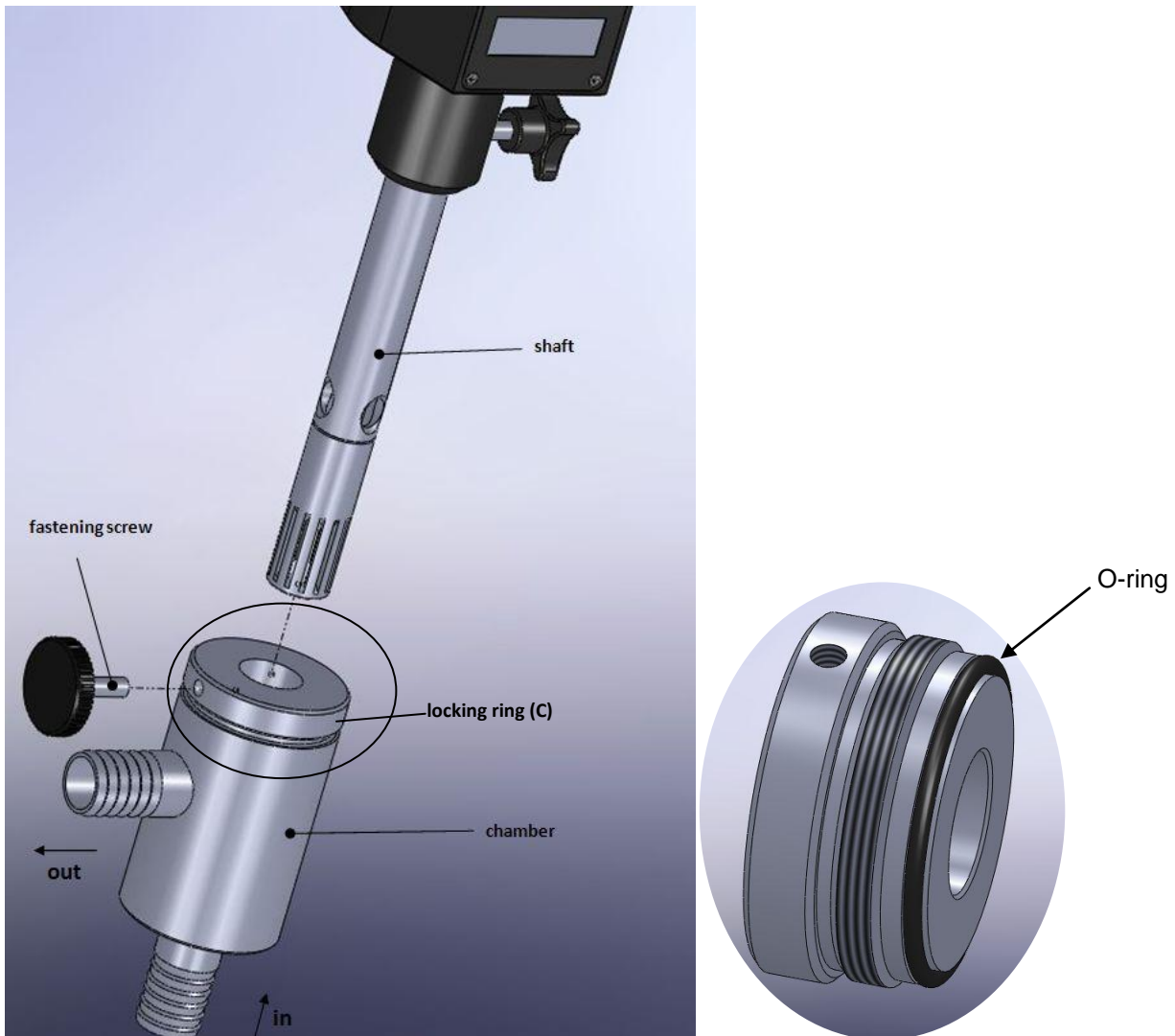Take care to avoid damage of parts during assembling respectively disassembling.

Only the Shafts G20 or G30 are suitable to be connected to the flow-through chamber DK 30; different Generator versions are available (F, N, V and M).

- Remove locking ring (C), now you will see a spacing washer and an O-ring
- Remove thrust washer and O-ring
- Now assemble O-ring, thrust washer and locking ring (C) onto Shaft G20 or G30
- Be sure O-ring and spacing washer are in place -
- Insert Generator end of Shaft G20 or G30 into chamber
- Carefully slide down O-ring, thrust washer and locking ring
- Screw locking ring in completely, to insure complete sealing -
- Push in Shaft G20 or G30 until it reaches the end of the chamber -
- Tighten fastening screw to secure Shaft G20 or G30. The inlet tube at the front end is connected to the system. From the outlet tube at the side the product is pumped into a container. It is possible to re-circulate, if necessary.

## 6 Operation of Flow Through-Chamber

### 6.1 Flow-Through Mode

Connect the vessel containing the media to the inlet by a tube (inner Diameter = 20 mm). Connect another tube (inner Diameter = 20 mm) to the outlet and lead it to a collecting basin.

**Note:** Secure all tube connections with suitable tube clips.

### 6.2 Circulation Mode

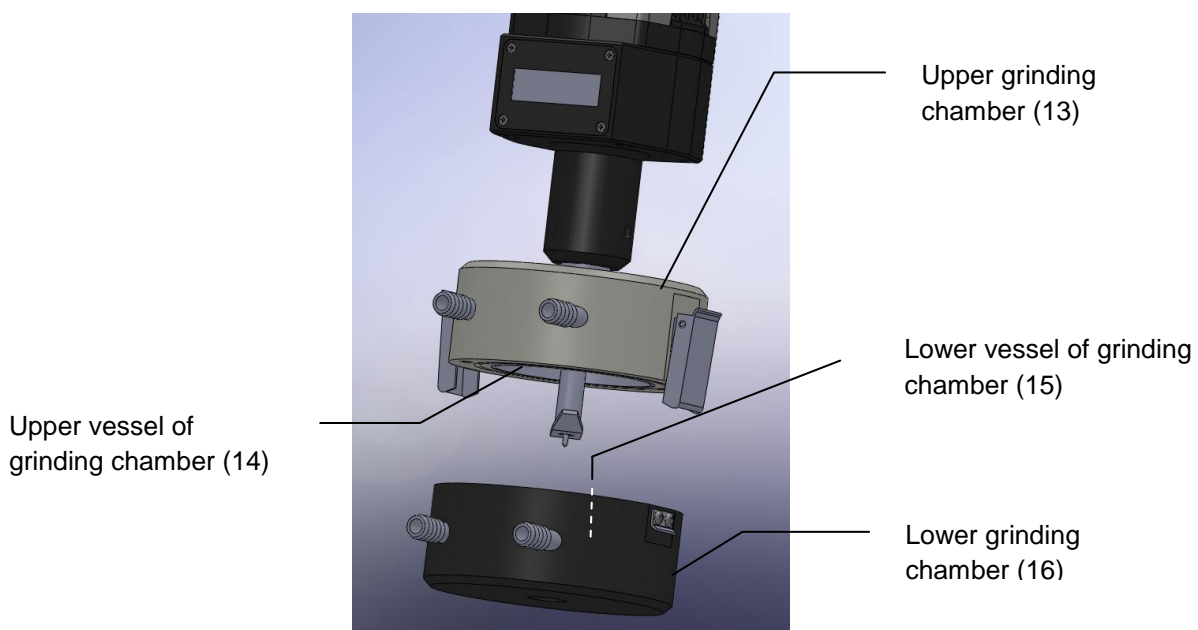
Connect in and outlet tube as described in 6.1 and lead the outlet tube back into the original vessel. This way the medium might flow several times through the in-line unit.

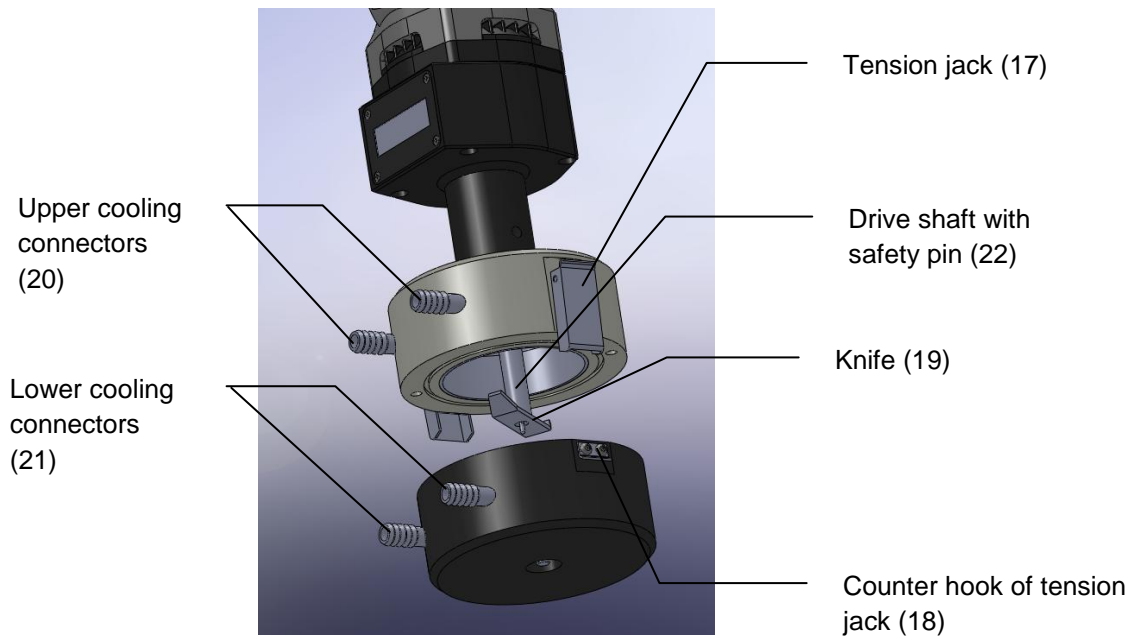
**Note:** The flow-through-chamber is not self-priming. Inlet tube and flow through chamber have always to be filled with a medium. We therefore recommend to set up the vessel containing the sample at the same height as the flow through chamber and to additionally fit a blockage (for example a ball valve) to the connecting tube.

## 7 Analytical Mill Attachment AX 60

### 7.1 Description of the Analytical Mill Attachment AX 60

For analysis, material testing and quality control of organic and inorganic substances of various hardness. The grinding chamber consists of two stainless steel beakers, both with separate cooling facilities. The total capacity of the grinding chamber is 180 cm<sup>3</sup>. The two grinding chambers are – when closed – sealed with an o-ring to prevent from dust and losses. This analytical mill is especially developed to also enable the use of coolants such as N<sub>2</sub> or to crush and analyse material that is best crushed when super cooled. The complete grinding chamber is connected by a mechanical clutch to the drive unit UNIDRIVE X1000 or UNIDRIVE X1000D. The speed control of these motors allow for a smooth start of the grinding mill at low speed. Extremely rapid disintegration of the material is possible due to the very high maximum speed of these motors. The grinding cutter knife (19) declutches when the grinding chamber is opened. The design prevents from unintended run of the cutter when the chamber is not closed.

## 8 Mounting of the Analytical Mill AX 60

- Make sure that the switch of the drive unit is in off position
- Close the chamber and connect it to the drive unit – therefore insert the drive shaft with the safety pin (22) into the coupling of the drive unit. Make sure that the upper part is properly fitted to the drive unit. This can be easily detected by the "click" that is heard when the coupling Shaft has reached the end of the coupling slot of the motor.
- Now fix the mill attachment firmly with the knurled screw at the coupling of the drive unit.
- Start the drive unit at the lowest speed.

**Important Note:** *There are 4 connections for the separate cooling facilities of the two grinding chambers. When using N<sub>2</sub> or other super cooling media please make sure that knife shaft and bearings do not freeze. This could occur due to condensed water.*

## 8.1 Exchange of the cutting knife

**Attention:**

*The cutting knife has sharp edges. Handle with care! Risk of injuries!*

- Prior to exchange the cutting knife disconnect the unit from the mains supply by unplugging the unit.
- Loosen the holding screw (6) and firmly pull out the mill attachment by hand. It will disconnect easily.
- For undoing the cutting knife (19), use the socket wrench (to be ordered separately). It is used to counter-hold the drive shaft (22) to undo the cutting knife.
- Turn the knife to the right
- Now screw on the exchange knife.
- Close the mill attachment and attach it to the motor coupling of the drive unit. Make sure that the coupling fits well. This can be easily detected by the "click" that is heard when the coupling shaft has reached the end of the coupling slot of the motor.
- Securely fasten the holding screw (6).

## 9 Maintenance and Cleaning

### 9.1 Cleaning the Drive Unit UNIDRIVE X1000/X1000D

To clean the drive unit use only water with a detergent that contains tensides or use isopropyl alcohol for stubborn soiling.

### 9.2 Cleaning the Homogenising Tool

*The Generators have sharp edges. Handle with care. Risk of injuries!*

To avoid clogging of the shaft, clean shaft, generator and in case of G-shaft the sealing parts after each use. This is done by operating it in a solvent which dissolves substance residues and is not harmful to the gasket. This is usually sufficient to clean the Generator.

**Chemical sterilization** may be also a method. General-purpose disinfectants such as formalin, alcohol, etc. may be used. It is important to remove disinfectant residues with sterilized water.

**Important Note:**

*Make sure that the bearings, O-rings and gaskets are resistant to solvents.*

### **Sterilization with moist heat:**

This denotes the use of a steam jet pressurized to 2 bars at 120°C.

### 9.3 Rough Cleaning of Flow-Through-Chamber and Tubing

Flow-through-chamber and tubing might be roughly cleaned by pumping a suitable liquid through the chamber instead of the medium.

### 9.4 Cleaning the Inside of the Flow-Through-Chamber

Before cleaning please disconnect first drive unit from mains and detach the flow through chamber DK 30

The flow-through chamber can be disassembled without disconnecting the tubes. **Attention:** Rest liquid escapes!!!

Please take care that neither O-rings nor any other parts experience any damage during assembling respectively disassembling the flow-through chamber.

**Attention:** Before operation check the slip ring sealing. The slip ring should be easily turned by hand. If it is stuck after a long period without use just loosen slip ring by wetting it with water. Never operate the flow-through chamber if the slip ring is stuck!!

**Note:** Rotor and stator wrenches are listed in the price list under accessories. Always unscrew the rotor before the stator.

### 9.5 Cleaning of the Analytical Mill Attachment AX 60

Before cleaning please disconnect first drive unit from mains and detach the mill attachment AX 60.

Please take care that neither the o-ring nor any other parts experience any damage during assembling respectively disassembling the mill attachment.

Remove residues from the inside of the chamber. To avoid clogging, clean shaft, knife and the inside of the chambers after each use. This is done by operating the attachment using a solvent which dissolves substance residues and is not harmful to the gasket. This is usually sufficient to clean the attachment inside.

**Chemical sterilization** may be also a method. General-purpose disinfectants such as formalin, alcohol, etc. may be used. It is important to remove disinfectant residues with sterilized water.

**Important Note:**

*Make sure that the bearings, O-rings and gaskets are resistant to solvents.*

**Sterilization with moist heat:**

This denotes the use of a steam jet pressurized to 2 bars at 120°C.

### 9.6 Maintenance of drive units UNIDRIVE X1000/X1000D

**Attention:**

*Do not open the instrument. Repairs are only to be carried out by trained service technicians.*

## 9.7 Maintenance homogenizing tools

The Generators have sharp edges. Handle with care. Risk of injuries!

The gaskets in the homogenizing tools must be constantly monitored. In the event of leakage the suction effect of the rotating Shaft can cause the medium to penetrate as far as the drive unit. If liquid emerges from the side hole at the top of the Shaft tube stop work immediately and check the gaskets. The function of the Generators depends on the condition of the sharp edges on the rotor and stator. These edges may be blunted very quickly in abrasive media, reducing the effectiveness of homogenizing.

## 10 Dismantling, Transport and Storage

### 10.1 Dismantling

1. Switch the unit off.
2. Disconnect it from the mains supply.
3. Remove any glass beakers and any other equipment around the drive unit.
4. Remove the homogenizing tool by loosening the holding screw (6).
5. Loosen the special clamp and remove the clamp from the support rod.
6. Unscrew the support rod.
7. Now you may remove the instrument from the working area.

### 10.2 Transport and Storage

#### Prior to transport:

Switch the instrument off and proceed with dismantling as described under “Dismantling”.

Place the instrument and its parts in its original packaging or another suitable container to protect it during transport. Close the packaging with adhesive tape.

Store the instrument in a dry environment. Please observe the specified conditions of the ambient (temperature and humidity).

Do not subject the instrument to mechanical shocks or vibration during transporting it.

In case you do not use the original packaging please mark the box with the following notes:

- Glass symbol (handle with care, fragile)
- Umbrella (keep dry)
- Content (list of content)
- Storage ambient:
  - Max. ambient temperature : RT to +40°C
  - Max. humidity: 80%

## 11 Disposal

Please dispose of used instruments and defective components at your local recycling collection point. Prior to disposal, sort according to materials: Metal, glass, plastic, etc. Also be sure to dispose of the packing material in an environmental-friendly manner.

## 12 Warranty and Liability

The manufacturer agrees to either repair, or replace, at the manufacturer's discretion, any defects in materials or workmanship which develop within 24 months of the delivery of this product to the original user. In the event of replacement, the replacement unit will be guaranteed for the remainder of the original twenty-four (24) months period or ninety (90) days, whichever is longer.

If this product should require service, contact your local distributor or manufacturer for necessary instructions.

This guarantee will not apply if the defect or malfunction was caused by accident, neglect, unreasonable use or fitness for a particular purpose, which extend beyond the description and period set forth herein.

The manufacturer's sole obligation under this guarantee is limited to the repair or replacement of a defective product and the manufacturer shall not, in any event, be liable for any incidental or consequential damages of any kind, resulting from use or possession of the product.

### **Attention:**

The user has to determine, if the instrument is suitable for his specific application. If there are any further queries, contact your local dealer or the manufacturer direct.

## 13 Technical Data

### 13.1 Technical Data UNIDRIVE X1000/X1000D

Type: UNIDRIVE X1000	Specifications
Motor Power intake	1050 Watt
Motor Power output	650 Watt
Idle speed	4000 - 33000 rpm
Voltage	230V/50 Hz or 110V/60 Hz
Case dimensions	75x75x290 mm, WxDxH
Speed Display	-
Weight	2.1 kg
Permissible ambient temperature	5 - 40 °C
Permissible humidity	80 % RH

Type: UNIDRIVE X1000D	Specifications
Motor Power intake	1050 Watt
Motor Power output	650 Watt
Idle speed	4000 - 33000 rpm
Voltage	230V/50 Hz or 110V/60 Hz
Case dimensions	78x75 x 310 mm, WxDxH
Speed Display	Yes
Weight	2.2 kg
Permissible ambient temperature	5 - 40 °C
Permissible humidity	80 % RH

## 14 Repairs

*When returning instruments for repair that have come into contact with hazardous substances, please:*

*Fill in attached "Repair Return Form"*

*Provide precise information on the relevant medium*

*Take protective measures to ensure the safety of our receiving and maintenance personnel*

*Mark the package as appropriate for hazardous materials.*

## 14.1 Repair Return Form

### CONTACT/USER INFORMATION

Contact:	Phone No.
Fax No.	Email:
Billing:	Shipping:
Company	Company
Address	Address

### INSTRUMENT INFORMATION

Model	Serial No.	
Please describe all problems/malfunctions		
Operating Conditions (please fill in if applicable)		
Ambient Temp.	Humidity	Speed
Load	Volume	Viscosity
Temperature in °C	Sample Temperature	Operating Time
Sample Description*		

\*NOTE: If the instrument was exposed to hazardous material, it must be decontaminated BEFORE returning it to CAT and an MSDS for hazardous material must be included with the instrument.

### RETURN SHIPPING

UPS <input type="checkbox"/>	Air Parcel Post <input type="checkbox"/>	Collect** <input type="checkbox"/>	Other** <input type="checkbox"/>
------------------------------	------------------------------------------	------------------------------------	----------------------------------

\*\*Your account number is required for UPS collect respectively the address and contact of your preferred forwarder if you choose any other transport means.

### PACKAGING INSTRUCTIONS TO RETURN A CAT INSTRUMENT FOR REPAIR

- ✓ Remove all accessories (e.g. homogniser tools, stirring paddles) from the instrument
- ✓ Clean excess testing material off the instrument/accessory
- ✓ Include MSDS sheets for all hazardous materials used with this instrument
- ✓ Pack the instrument in its original box. If the box is not available, take care to wrap the instrument and accessories with enough material to support them.
- ✓ DO NOT send pedestal stand unless there is a problem with the upright rod, clamp or base. If there is a problem with the stand remove the upright rod from the base and individually wrap each item to avoid contact with the instrument. (Applicable for overhead stirrers and homogenisers.)
- ✓ Pack the instrument and related items in a strong box for shipping. Mark the outside of the box with handling instructions.

Example: "Handle with care" or "Fragile- Delicate Instrument"

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